


HARDNESS TEST

Article Nr. T1454-00

TEST PROCEDURE

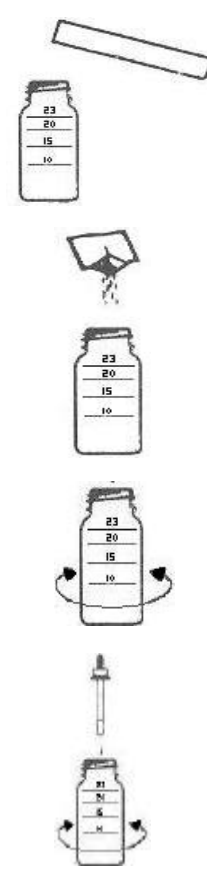

1. Rinse plastic tube with sample water before test. Fill a plastic tube full with sample water.
Pour the volume of the plastic tube in the mixing bottle.
2. Add the content of one Uni Ver[®] 3 Hardness Reagent Powder Pillow to the mixing bottle.
3. Swirl to mix.
4. While swirling the mixing bottle, add Hardness 3 Titrant solution drop by drop.
5. When the sample changes from red to blue, record the number of drops added. The hardness equals the number of drops of Hardness 3 Titrant solution added in.

Drops = Hardness



MARITECH- MARINE TECHNOLOGIES AB

P.O.Box 143, SE-291 22 Kristianstad, SWEDEN · Tel. Int. +46 44 106040, Fax. Int. +46 44 106440
e-mail: marisol@maritech.se www.marisol.nu